



PID Orientation Tube

Ref: CYT-PIDOT

RUO

For Research Use Only. Not for use in diagnostics procedures

VIALS OF PID Orientation Tube KIT ARE A LYOPHILIZED PRODUCT. READ CAREFULLY THE FOLLOWING INSTRUCTIONS FOR RECONSTITUTION:

The PID Orientation Tube preserves the stability of the pre-mixed combination of antibodies until expiration date shown on the label. Reconstitute each lyophilized vial containing the pre-mixed combinations with 300 µl of distilled water. Unused volume of reconstituted vial is stable during four weeks from date of reconstitution if it is stored in the dark at 2-8° C.

INTENDED USE

PID Orientation Tube kit contains 4 lyophilized vials. Each vial contains a 10 conjugated antibodies mixture designed for the screening of primary immunodeficiencies by flow cytometry (FC). This 6-colors tube can be used for the study of defects and numerical changes in the relative distribution of the different leukocyte populations and subpopulations ⁽¹⁻³⁾ (antigen presenting cells, memory, naïve and effector B and T cells, etc.). This reagent must be used by FC qualified personal.

REAGENT COMPOSITION

Material included

PID Orientation Tube kit contains sufficient volume for 20 tests and includes:

- Following reagents for specific tandem compensation ready to use in liquid format sufficient for 5 test (5µl/test):
 - CD45-APC-C750. Specific compensation reagent for APC-C750 fluorochrome.
 - CD4-PerCP-Cyanine5.5 (Specific compensation label CD4/IgM-PerCP-Cyanine5.5). Specific compensation reagent for PerCP-Cyanine5.5.
 - CD19-PECyanine7 (Specific compensation label CD19/TCRγδ-PE-Cyanine7). Specific compensation reagent for PECyanine7.
- A combination of Surface Membrane (Sm) staining antibodies distributed in 4 lyophilized vials of 5 tests each and identified as follows:
 - Antibody anti human IgD/CD8-FITC, clone: IADB6/UCHT-4, isotype: IgG2a/IgG2a
 - Antibody anti human CD16/CD56-PE, clone: 3G8/C5.9, isotype: IgG1/IgG2b
 - Antibody anti human CD4/IgM-PerCP-Cyanine5.5, clone: RPA-T4/MHM-88, isotype: IgG1 /IgG1
 - Antibody anti human CD19/TCRγδ-PE-Cyanine7, clone: 19-1/TCR-1, isotype: IgG1/IgG1
 - Antibody anti human CD3-APC, clone: UCHT-1, isotype: IgG1
 - Antibody anti human CD45-APC-C750, clone: HI30, isotype: IgG1

Fluorochrome	FITC		PE		PerCP-Cyanine5.5		PE-Cyanine7		APC	APC-C750
	IgD	CD8	CD16	CD56	CD4	IgM	CD19	TCR-γδ	CD3	CD45
Marker	IgD	CD8	CD16	CD56	CD4	IgM	CD19	TCR-γδ	CD3	CD45
Clone	IADB6	UCHT-4	3G8	C5.9	RPA-T4	MHM-88	19-1	TCR-1	UCHT-1	HI30
Isotype	IgG2a	IgG2a	IgG1	IgG2b	IgG1	IgG1	IgG1	IgG1	IgG1	IgG1

- Fixative free ammonium chloride erythrocyte lysing solution (BulkLysis™).

Material required but not included

- Reagents for generic compensation of non-tandems (Cytognos recommends to follow the EuroFlow instrument set-up and compensation ⁽⁴⁾).
- Rainbow beads calibration particles ⁽⁴⁾.
- Antibodies against human CD27-BV421 and CD45RA-BV510 that must be used as drop in.
- Test tubes suitable for obtaining samples in the flow cytometer used. Usually tubes with a rounded bottom for 6 ml, 12x75 mm are used (FACS tubes).
- 50 ml tubes (Falcon).
- Automatic pipette and tips.
- Micropipette and tips.
- Vortex mixer.
- Roller mixer.
- Chronometer.

- Centrifuge.
- Pasteur pipette or vacuum system.
- Distilled water.
- Erythrocyte fixative lysing solution.
- Wash buffer as phosphate buffered saline (PBS) + 0,09% (m/v) of NaN₃ + 0,5% (m/v) of Bovine Serum Albumin (BSA).

STORAGE CONDITIONS

PID Orientation Tube kit is stable until the expiration date shown on the label, when stored at 2-8° C. The expiration date applies to the lyophilized product. After reconstitution vials with the pre-mixed combination are stable for four weeks when stored at 2-8° protected from light.

Components should not be frozen or exposed to direct light during storage or during incubation with cells. Keep vials in a dry place. Once opened, vials must be stored in a vertical position to avoid any possible spillage.

WARNINGS AND RECOMMENDATIONS

1. For research use only. Not for use in diagnostic procedures.
2. If components of this kit are altered by addition of other components, such conditions must be validated by the user.
3. The kit is stable until the expiration date shown on the label if it is properly stored. Do not use after the expiration date shown on the label. If the reagents are stored in conditions different from those recommended, such conditions must be validated by the user.
4. Alteration in the appearance of the reagents, such as the precipitation or discoloration indicates instability or deterioration. In such cases, the reagents should not be used.
5. It contains 0,09% (m/v) sodium azide (CAS-No. 26628-22-8) as a preservative, but even so care should be taken to avoid microbial contamination of reagent or incorrect results may occur.

Indication(s) of danger:

H302 Harmful if swallowed

Safety advice:

P264 Wash thoroughly after handling.

P270 Do not eat, drink or smoke when using this product.

P301+P312 If swallowed, call a poison center or doctor/physician if you feel unwell.

P301+P330 If swallowed, rinse mouth.

P501 Dispose of contents/container in accordance with local/regional/national/international regulation.

6. All patient specimens and materials with which they come into contact are considered biohazards and should be handled as if capable of transmitting infection⁽⁶⁾, and disposed according to the legal precautions established for this type of product. Also recommended is handling of the product with appropriate protective gloves and clothing, and its use by personnel sufficiently qualified for the procedures described. Avoid contact of samples with skin and mucous membranes. After contact with skin, wash immediately with plenty of water.
7. Use of the reagents with incubation times or temperatures different from those recommended may cause erroneous results. Any such changes must be validated by the user.
8. Any serious incident relating to the product must be reported to Cytognos SL as well as the competent professional authority of the Member State in which the user is established.

PROCEDURE

Reconstitute each lyophilized vial containing the pre-mixed combinations with 300µl of distilled water, mix well and let the solution at least 30 minutes before using.

Preparation

Since some cellular subpopulations have a very low frequency in peripheral blood, it is recommended to use the macro-volume lysis technique with a hypotonic solution without fixative, prior to labeling.

Sample must be collected in commercially available anticoagulant-treated tube (use of EDTA is recommended)^(6, 7).

1. Determine the absolute count of leukocytes per µl for the sample to be processed.
2. Transfer 1.5mL (nucleated cells > 8x10⁶/ml) or 2.0ml (nucleated cells < 8x10⁶/ml) of sample in a 50 ml-Falcon tube.
3. Do not use more than 2 ml of sample per 50 ml of lysing solution. If larger volumes of sample need to be processed (i.e. starting cells concentration is low), use several 50 ml tubes.
4. Fill the tube up to reach 50 ml volume with BulkLysis™ (CYT- BL) (diluted to 1X in distilled water and at room temperature - RT).
5. Mix well and incubate for 15 min in a roller or sample-shaker device.
6. Centrifuge at 800 g for 10 min and remove the supernatant using a Pasteur pipette or a vacuum system without disturbing the cell pellet. Typically, 300 µl of cell suspension should remain in the tube.
7. Add 2 ml of PBS + 0,09% (m/v) of NaN₃ + 0,5 % (m/v) of BSA and resuspend the cell pellet vigorously.
8. Complete the volume of the tube containing the cell suspension up to 50 ml final volume with PBS + 0,09% (m/v) of NaN₃ + 0,5 % (m/v) of BSA.
9. Mix well.
10. Centrifuge at 800 g for 5 min and remove the supernatant using a Pasteur pipette or a vacuum system, without disturbing the cell pellet.
11. Resuspend the cell pellet in 2 ml of PBS + 0,09% (m/v) of NaN₃ + 0,5 % (m/v) of BSA. Mix well and transfer this volume to a 5 ml "FACS tube".

12. Wash the 50 ml Falcon tube with 2 mL of PBS + 0,09% (m/v) of NaN₃ + 0,5 % (m/v) of BSA more to recover cells that might have left in the original tube. Add this volume to the 5 ml tube containing the rest of the sample transferred in step 10.
13. Centrifuge at 540 g for 5 min and remove the supernatant by decanting or using a Pasteur pipette. If the remaining cell volume is lower than 300 µl, PBS + 0,09% (m/v) of NaN₃ + 0,5 % (m/v) of BSA will be added to reach a volume of at least 300 µl.
14. In case multiple 50 ml tubes were used (because it was needed to lyse large sample volumes) the cell suspensions from the same sample should be combined at this moment, before adjusting cell concentration. Try to keep the final volume low, so that, in case that cell concentration needs to be adjusted as indicated in the next step, it can be easily done by diluting with the recommended buffer.
15. Adjust the final cells concentration to 1×10^5 cells/ µl, by resuspending the pellet with PBS + 0,09% (m/v) of NaN₃ + 0,5 % (m/v) of BSA.
16. Adjust the volume in order to obtain 50 µl containing 2.5×10^6 cells of the cell suspension per each tube to be stained/acquired. Following this procedure there will be an excess of sample that could be used in other panels, according to the purpose.

Surface Staining:

1. Add 50 µl of antibody mixture in the tube containing the 50 µl cell suspension (2.5×10^6 cells per tube). Add proper volume of CD27 and CD45RA (not included in the mixture).
2. Mix well. For optimal staining conditions, complete with PBS until a final volume of 200 µl.
3. Incubate for 30 min at RT protected from light.
4. Add 2 ml of fixative Lysing Solution.
5. Mix well.
6. Incubate for 10 min at RT protected from light.
7. Centrifuge for 5 min at 540 g.
8. Discard the supernatant using a Pasteur pipette or vaccum system without disturbing the cell pellet, leaving approximately 100 µl residual volume in each tube.
9. Add 2 ml of PBS+0,5% (m/v) BSA + 0,09% (m/v) NaN₃ to the cell pellet.
10. Mix well.
11. Centrifuge for 5 min at 540 g.
12. Discard the supernatant using a Pasteur pipette or vaccum system without disturbing the cell pellet, leaving approximately 100 µl residual volume in each tube.
13. Resuspend the cell pellet in 300 µl PBS + 0,5% (m/v) BSA (without NaN₃).
14. Acquire the cells immediately after staining or, if not immediately acquired, store at 4°C, for 1h maximum, until measured in the flow cytometer.
15. Acquire the sample at medium flow rate (at least 1×10^6 cells).

LIMITATIONS

- Blood samples should be stored at 18-22°C and be tested within the 24 hours after they are obtained.
- It is advisable to acquire stained samples on the cytometer as soon as possible to optimize the results. Non viable cells may stain nonspecifically. Prolonged exposure of whole blood samples to lytic reagents may cause white cell destruction and loss of cells from the target population.
- When using whole blood procedures, all red blood cells may not lyse under following conditions: nucleated red blood cells, abnormal protein concentration or hemoglobinopathies. This may cause falsely decreased results due to unlysed red blood cells being counted as leucocytes.
- Results obtained by FC may be erroneous if the cytometer laser is misaligned or the gates are improperly set.
- Each laboratory should establish a normal range for lymphocyte subsets. We recommend to follow the EuroFlow antibody panels together with the EuroFlow instrument set-up, sample preparation and data analysis procedures ⁽⁴⁾.
- Cells separated from whole blood by means of density gradients may not have the same relative concentrations of cells as unseparated blood. This may be relatively insignificant for samples from individuals with normal white blood cell counts. In leucopenic patients, the selective loss of specific subsets may affect the accuracy of the determination.

QUALITY CONTROL

- Pipettes precision and cytometer calibration should be verified to obtain optimal results.
- In multicolour panels, fluorochromes emit in wavelengths that can show certain spectral overlap which must be corrected by electronic compensation. Optimal compensation levels can be established by analysing cells from healthy individuals stained with mutually exclusive monoclonal antibodies conjugated with appropriate fluorochromes.
- This product has been manufactured in accordance with standards of production and quality system of the ISO 9001:2008 standard.

REFERENCES

1. Sigal LH. Basic science for the clinician 58: IgG subclasses. J Clin Rheumatol. 2012 Sep;18(6):316-8.
2. Warnatz K1, Denz A, Dräger R, Braun M, Groth C, Wolff-Vorbeck G, Eibel H, Schlesier M, Peter HH. Severe deficiency of switched memory B cells (CD27(+)/IgM(-)/IgD(-)) in subgroups of patients with common variable immunodeficiency: a new approach to classify a heterogeneous disease. Blood. 2002 Mar 1;99(5):1544-51.
3. The Immune Response. Basic and Clinical Principles (2006). Tak W. Mak and Mary E. Saunders. Elsevier Academic Press.
4. Kalina T, Flores-Montero J, van der Velden VHJ, Martin-Ayuso M, Böttcher S, Ritgen M, et al. EuroFlow standardization of flow cytometry instrument settings and immunophenotyping protocols. *Leukemia*, 26(9):1986-2010 (2012).

5. Protection of Laboratory Workers from occupationally acquired infections. Second edition; approved guideline (2001). Villanova PA: National Committee for Clinical Laboratory Standards; Document M29-A2.
6. Procedures for the collection of diagnostic blood specimens by venipuncture- approved standard; Fifth edition (2003). Wayne PA: National Committee for Clinical Laboratory Standards; Document H3-A5.
7. Clinical applications of flow cytometry: Quality assurance and immunophenotyping of lymphocytes; approved guideline (1998). Wayne PA: National Committee for Clinical Laboratory Standards; Document H42-A.

WARRANTY

This product is warranted only to conform to the quantity and contents stated on the label. There are no warranties that extend beyond the description on the label of the product. Cytognos's sole liability is limited to either replacement of the product or refund of the purchase price.

EXPLANATION OF SYMBOLS

	Use by (YYYY-MM)
	Storage temperature limitation
	Keep out of sunlight
	Consult instructions for use
	For research use only
	Batch code
	Catalogue number
	Manufacturer

PRODUCED BY **CYTOGNOS SL**

Polígono La Serna, Nave 9
37900 Santa Marta de Tormes
Salamanca (España)

Phone: + 34-923-125067

Fax: + 34-923-125128

Ordering information: admin@cytognos.com

Technical information: support@cytognos.com

www.cytognos.com